

uvex

Disposable Protective Clothing

protecting people










uvex

uvex 5/6 air



Disposable Protection Clothing


Guide to standards and products

	uvex 5/6 air	uvex 5/6 air	uvex 5/6 comfort light	uvex 5/6 comfort	uvex 5/6 classic light	uvex 5/6 classic	uvex 4B	uvex 3B chem light	uvex 3B chem classic	uvex 3B chem classic*
	98173	89976	88497	98710	17595	98449	98375	89843	89880	88498
EN 14605  Liquid-tight coveralls								■	■	■
EN 14605  Spray-tight coveralls							■	■	■	■
EN 13034  Limited spray-tight coveralls (protection against liquid mist)	■	■	■	■	■	■	■	■		
ISO 13982-1  Particle-tight coveralls (protection against solid particles)	■	■	■	■	■	■	■	■	■	■
EN 1149-5  Protective clothing with antistatic properties (antistatic function only guaranteed if air humidity > 25%)	■	■	■	■	■	■	■	■	■	■
EN 14126  Protective clothing against infective agents					■	■	■	■	■	■
EN 1073-2  Protective clothing against particulate radioactive contamination (no protection against radioactive rays)	■	■	■		■	■	■	■	■	■
DIN 32781  Protective clothing against pesticides							■			
ISO 27065  Protective clothing for use with pesticides								■		

For more detailed information, please see the relevant certificates, technical data sheets or the endorsement on the product!

Defining environmental conditions

Are you exposed to dusty, damp or wet conditions during your work? To make it easier for you to select the right clothing for the job, we've categorised our overalls depending on their suitability for these environments. The bars indicate how suitable the product is for each setting. We also provide information on breathability so that you can take wearer comfort into account in your decision.

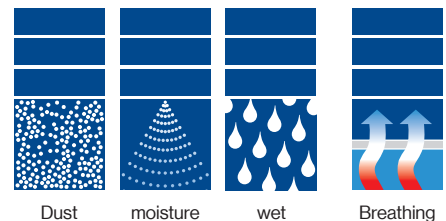

Identifying additional characteristics

In some work environments, there are additional requirements that you will need to take into account. These symbols indicate whether the items are free from silicone and paint wetting impairment substances.



Disposable protective clothing

Type 4 and Type 5/6 features and technologies

1

3 pieced hood

- optimum fit
- no limitations of view
- perfect match with respirators and goggles
- soft elastic on edge, feels comfortable

2

Zipper flap

- Zipper flap with integrated self adhesive tape
- easy to gum up
- additional security

3

Fingerloop

- finger loops made of soft elastic
- to keep the sleeves at the wrist
- additional security

4

Elastic in waistband

- integrated elastic in waist for optimal fit
- ideal wearer comfort

5

elasticated bands on arm and legs

- convenient closures thanks to elastics at wrist and ankles for additional security
- soft elastic on edge for high wearer comfort



Technologies




uvex SMS

- build out of several layers of polypropylene non woven material
- Mixture of spunbond and melt-blown non woven material for a perfect balance between stability and softness
- High degree on air permeability and breathability



uvex com4-Laminat

- non woven polypropylene material laminated with a microporous polyethylene film
- non woven material on the inside for a comfortable feeling
- material is not sticking on skin
- the microporous film provides a certain level on breathability

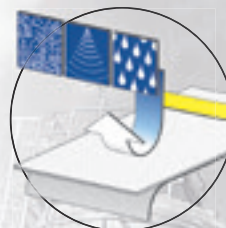

uvex Laminat light

- non woven polypropylene material laminated with a polyethylene film
- non woven material on the inside for a comfortable feeling
- material is not sticking on skin
- using the film is providing a higher level of protection, but reduces the breathability



uvex Laminat classic

- non woven polypropylene material laminated with a polypropylene film
- non woven material on the inside for a comfortable feeling
- material is not sticking on skin
- the polypropylene film provides an effective barrier against many organic and inorganic liquid chemicals

uvex 3B chem classic

Disposable coverall chemical protection Type 3B



89880



88498

Available from
January 2022

uvex 3B chem classic • uvex 3B chem classic+

General features:

- highly durable material combined with ultrasonically welded and taped seams ensure a great barrier and maximum of safety
- lightweight and durable material
- skin-friendly textile grip on inside

Protection features:

- offers protection against a wide range of chemicals
- optimal protection thanks to self-adhesive zip-per flap

Comfort features:

- elasticated waistband, elasticated bands on hood, arms and legs for a perfect fit
- middle finger loops prevent sleeves sliding up the arm

Areas of application:

- handling organic and inorganic chemicals
- cleaning and maintenance work
- chemical and pharmaceutical industries
- food industry
- remediation of soil contamination and dismantling
- industrial cleaning and maintenance
- tank cleaning
- work with paints and varnishes
- disposal of hazardous materials
- agricultural industry
- waste water treatment and drainage construction
- waste management
- disaster response and emergency services
- veterinary medicine and disease control
- oil and petrochemicals

Details for 88498:



The double zip flap makes it easy to put on



Double sleeves for extra safety



integrated boot sock for increased safety

Certified in acc. with



EN 14126



EN 14605



EN 14405



ISO 13982-1

TYPE 3B



EN 1073-2




TYPE 4B

TYPE 5B



Type 3

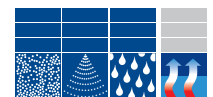

uvex 3B chem classic		
Art. no.	89880	88498
Material	polypropylene spunbond, laminated with polypropylene film	polypropylene spunbond, laminated with polypropylene film
Colour	yellow	yellow
Sizes	S to 3XL	S to 3XL
Order unit	PC	PC
Outer packaging	25 PC per CT	25 PC per CT



uvex-safety.com/en/overalls

uvex 3B chem light

Disposable coverall chemical protection Type 3B



89843

uvex 3B chem light

General features:

- light and highly flexible material for high wearer comfort with skin-friendly non-woven material inside

Protection features:

- ideal for cleaning work
- optimal protection thanks to self-adhesive zipper flap

Comfort features:

- elasticated waistband, elasticated bands on hood, arms and legs for optimal fit
- middle finger loops for securing the sleeves


Areas of application:

- handling low-concentration chemicals
- industrial and building cleaning
- ship building and automobile manufacture
- chemical and pharmaceutical industries
- handling paints and varnishes
- electronics
- handling and dismantling of asbestos
- remediation of contaminated sites
- livestock breeding and veterinary medicine
- waste management



Stitched and taped seams. Optimum protection against penetration by liquid aerosols and solid particles.

Certified in acc. with

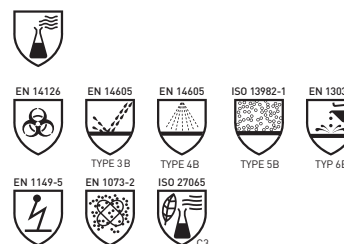

uvex 3B chem light	
Art. no.	89843
Material	polypropylene spunbond laminated with polyethylene film
Colour	white, yellow
Sizes	S to 3XL
Order unit	PC
Outer packaging	40 PC per CT

The chemicals included in this list are examples only. For a full overview for each individual garment, please consult the technical datasheets. You can also search our Chemical Expert System (see page 177).

Chemical	CAS no.	89843	89880
		Class in acc. EN 14325	Class in acc. EN 14325
Acetone	67-64-1	6 of 6	1 of 6
Acetonitrile	75-05-8	6 of 6	0 of 6
Iron(III) chloride (aq., 45%)	7705-08-0	–	6 of 6
Hydrofluoric acid (aq., 49%)	7664-39-3	6 of 6	6 of 6
Hydrofluoric acid (aq., 71 to 75%)	7664-39-3	–	5 of 6
Hydrofluoric acid in urea (62 to 64%)	7664-39-3	–	2 of 6
Formaldehyde (aq., 10%)	50-00-0	–	6 of 6
Isopropanol	67-63-0	6 of 6	6 of 6
Methanol	67-56-1	6 of 6	6 of 6
Sodium hydroxide, 50%	1310-73-2	6 of 6	6 of 6
Nitric acid, 69.5%	7697-37-2	–	6 of 6
Hydrochloric acid (aq., 37%)	7647-01-0	5 of 6	6 of 6
Sulphuric acid, 96%	7664-93-9	6 of 6	6 of 6
Hydrogen peroxide	7722-84-1	6 of 6	6 of 6

The information provided in the table was obtained in tests conducted under laboratory conditions (at temperatures of 21 ± 2°C). As additional stresses, such as higher temperatures and mechanical influences, are often present in practice, these values serve as a guide only. The information is provided without obligation and is not intended to replace your own suitability tests.

uvex

uvex 4B · uvex 3B chem light

Protection against pesticides – DIN 32781 / ISO 27065

In farming and forestry applications, workers are repeatedly exposed to hazardous chemical and biological substances.

Depending on the type of application, the duration of the exposure and the specific effects of the hazardous substance, it may be necessary to wear personal protective equipment.

In addition to respiratory protection and gloves, appropriate chemical protection clothing is an important component of the wearer's protective equipment.



You can find detailed information about this topic in our related blog post

uvex 4B

Disposable coverall chemical protection type 4B



98375



89094



Stitched and taped seams. Optimum protection against penetration by liquid aerosols and solid particles.



uvex 4B

General features:

- light and highly flexible material for high wearer comfort lined with skin-friendly non-woven material inside

Protection features:

- taped seams for optimal protection
- self-adhesive zipper flap for optimal protection


Comfort features:

- elasticated waistband, elasticated bands on hood, arms and legs for optimal fit
- middle finger loops prevent sleeves sliding up the arm

Areas of application:

- low-pressure industrial cleaning and building cleaning
- ship building and automobile manufacture
- chemical and pharmaceutical industries
- handling paints and varnishes
- agriculture and horticulture
- pest control
- electronics
- work with asbestos and dismantling
- remediation of contaminated sites
- pharmaceutical industry and laboratories
- sampling
- livestock breeding and veterinary medicine
- waste management

Certified in acc. with

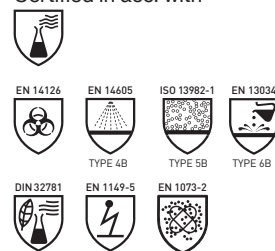

	uvex 4B	uvex 4B with overshoes
Art. no.	98375	89094
Material	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polyethylene film
Colour	white, orange	white, orange
Sizes	S to 3XL	S to 3XL
Order unit	PC	PC
Outer packaging	40 PC per CT	40 PC per CT



uvex-safety.com/en/overalls

uvex 5/6 classic

Disposable coverall chemical protection Type 5/6



98449



17595



uvex 5/6 classic

General features:

- extremely lightweight and breathable material guarantees an excellent combination of comfort and safety
- skin-friendly and soft non-woven material on the inside

Protection features:

- self-adhesive zipper flap for optimal protection

Comfort features:


- elasticated waistband, elasticated bands on hood, arms and legs for a perfect fit
- middle finger loops prevent sleeves sliding up the arm

Areas of application:

- handling chemicals in dust and powder form
- varnishing and protection against paint splashes
- fibreglass production and processing
- industrial cleaning and maintenance
- automotive industry
- grinding and polishing
- cement manufacture
- mining and quarrying
- demolition and refurbishment
- working with asbestos
- wood and metal processing
- construction industry
- pharmaceutical industry

► other articles on request

Certified in acc. with

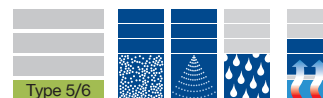

	uvex 5/6 classic	uvex 5/6 classic light
Art. no.	98449	17595
Material	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polyethylene film
Colour	white	white
Sizes	S to 3XL	S to 3XL
Order unit	PC	PC
Outer packaging	50 PC per CT	50 PC per CT



uvex-safety.com/en/overalls

uvex 5/6 comfort

Disposable coverall chemical protection Type 5/6



98710



88497

Available from
January 2022



98710: Breathable polyethylene laminate combined with an air permeable SMS back section to provide a high level of moisture management and breathability.



88497: Breathable polyethylene laminate combined with an air permeable SMS back section to provide a high level of moisture management and breathability.



uvex 5/6 comfort

General features:

- the combination of polyethylene laminate and a SMS back section provides a high level of moisture management, without reducing protection

Protection features:

- 98710: bound seams in contrasting colour
- 88497: with inside overlock seam
- self-adhesive zipper flap for optimal protection


Comfort features:

- optimal fit with elasticated waistband, elasticated bands on the hood, arms and legs
- middle finger loops prevent sleeves sliding up the arm

Areas of application:

- handling chemicals in dust and powder form
- varnishing and protection against paint splashes
- fibreglass production and processing
- industrial cleaning and maintenance
- automotive industry
- grinding and polishing
- cement manufacture
- mining and quarrying
- demolition and refurbishment
- working with asbestos
- wood and metal processing
- construction industry
- cleanroom applications
- pharmaceutical industry

Certified in acc. with

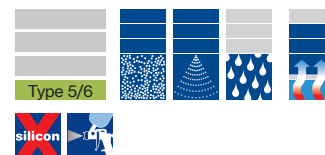

	uvex 5/6 comfort	uvex 5/6 comfort light
Art. no.	98710	88497
Material	Main material: polypropylene spunbond, laminated with polyethylene film back: SMS spunbond	Main material: polypropylene spunbond, laminated with polyethylene film back: SMS spunbond
Colour	white, lime	white
Sizes	S to 3XL	S to 3XL
Order unit	PC	PC
Outer packaging	40 PC per CT	40 PC per CT

uvex 5/6 air

Disposable coverall chemical protection Type 5/6



98173



89976



uvex 5/6 air

General features:

- excellent moisture management
- ideal for warm working environments and extended periods of wear

Protection features:

- optimal protection in dusty environments
- self-adhesive zipper flap for optimal protection

Comfort features:

- optimal fit with elasticated waistband, elasticated bands on hood, arms and legs
- middle finger loops prevent sleeves sliding up the arm

Areas of application:

- handling chemicals in dust and powder form
- working with asbestos
- fibreglass, ceramic fibre and synthetic resin production and processing
- wood and metal processing
- construction industry
- automotive industry
- grinding and polishing
- cement manufacture
- mining and quarrying
- refurbishment and renovation
- pharmaceutical industry
- repair work

Certified in acc. with



ISO 13982-1

EN 13034




TYPE 5

TYPE 6



EN 1149-5

EN 1073-2

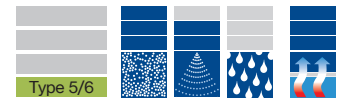

Type 5/6

uvex 5/6 air	
Art. no.	98173
Material	polypropylene spunbond
Colour	white
Sizes	S to 3XL
Order unit	PC
Outer packaging	50 PC per CT



uvex-safety.com/en/overalls

Disposable chemical protection

Accessories



9875200



9875300



9874946
9874948



9875046



8935346



8909746



8909546



8909646




Hood · Gauntlets · Overshoes · Overboots

	Hood	Gauntlets	Overshoes	Overboot
Art. no.	9875200	9875300	9874946 9874948	9875046
Sizes	one size fits all	one size fits all	42 to 46 46 to 48	42 to 46
Version	with velcro fastening	elasticated band on both ends	opening with elasticated band	opening with elasticated band and tie
Material	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polypropylene film
Colour	white	white	white	white
Order unit	BG	BG	BG	BG
Content	50 PC, loose	50 PR, loose	100 PR, loose	50 PR, loose

	Overboot	Overshoe with anti-slip	Overboots with anti-slip	Overboots with anti-slip
Art. no.	8935346	8909746	8909546	8909646
Sizes	42 to 46	42 to 46	42 to 46	42 to 46
Version	opening with elasticated band and tie	with anti-slip sole, elasticated opening improved fit	with anti-slip sole, elasticated opening with band improved fit	with anti-slip sole, elasticated opening with band
Material	polypropylene spunbond, laminated with polypropylene film	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polyethylene film	polypropylene spunbond, laminated with polyethylene film
Colour	yellow	white	white	yellow
Order unit	BG	BG	BG	BG
Content	30 PR, loose	100 PR, loose	25 PR, loose	30 PR, loose

Disposable chemical protection

Protection against infectious diseases

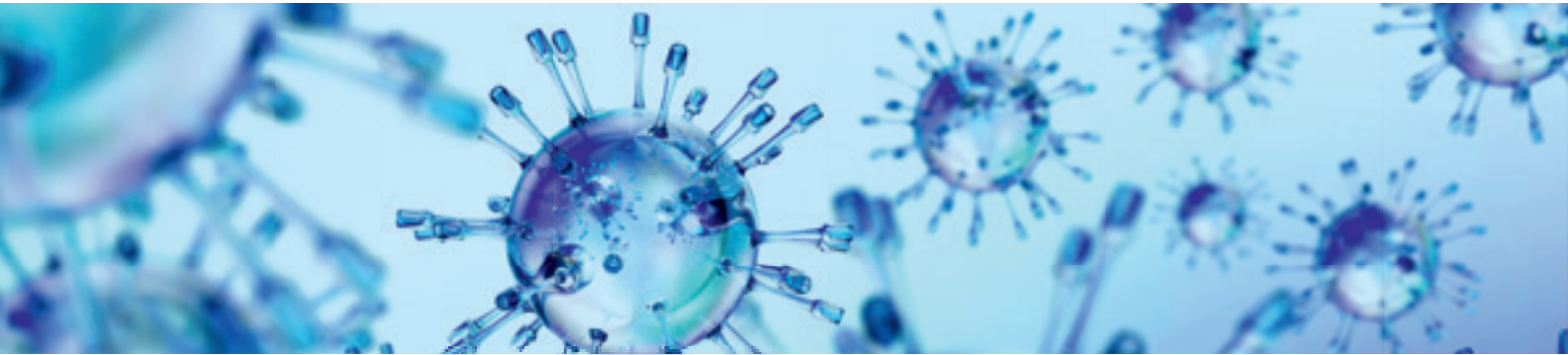

Workers may come into contact with biological substances in a wide range of situations. Appropriate personal protective equipment must be used to prevent potential infection.

Clothing designed to provide protection against infectious diseases establishes a physical barrier between the wearer's skin and the source of the infection, and prevents the disease from spreading to other people or environments.

Generally, micro-organisms such as bacteria, fungi and viruses are classed as biological substances. A detailed definition is provided in EU Directive 2000/54/EC. The crucial common denominator between all of these materials is that they can cause infections, trigger allergies, or produce a toxic effect.

Biological substances are placed in one of four categories depending on the infection risk they present:

Risk group 1:
Biological substances that are unlikely to cause disease in humans.

Risk group 2:
Biological substances that can cause human disease and might be a hazard to workers, but are unlikely to spread to the community; there is usually effective prophylaxis or treatment available.

Risk group 3:
Biological substances that can cause severe human disease and present a serious hazard to workers; they may present a risk of spreading to the community, but there is usually effective prophylaxis or treatment available.

Risk group 4:
Biological substances that cause severe human disease and are a serious hazard to workers; they may present a high risk of spreading to the community, and there is usually no effective prophylaxis or treatment available.

The risk groups of each biological substance are described in detail in the annex to EU Directive 2000/54/EC.

- Typical applications in which workers may be exposed to infectious substances include:
- Waste water treatment and drainage work
 - Waste disposal
 - Agriculture
 - Food industry
 - Work that involves contact with animals and/or products of animal origin
 - Healthcare, hospitals, emergency services

A detailed list of all biological substances that you may be exposed to in various applications, and the potential diseases associated with these substances, can be found in a number of publications, including the BGIA Report 1/2013.

Protective clothing in accordance with EN 14126:2003

The EN 14126 standard defines the performance requirements for clothing designed to protect against infectious diseases. The defined test procedures focus on the medium that contains the micro-organisms, such as fluids, aerosols or solid dust particles. Due to the heterogeneous nature of the micro-organisms, it is not possible to define performance criteria for them.

The tests stipulated by the standard relate only to the material; seam technologies are not taken into account. Taped seams provide a higher level of protection, as the micro-organisms are small enough to penetrate through the tiny needle holes along the seam. The entire protective suit must also be certified as chemical protection clothing (see page 177).

Wearing appropriate protective clothing does not provide complete, guaranteed protection against all chemical risks. It is also essential that you put on and remove the PPE correctly to ensure safety (see page 178). Any person who assists in the removal of the clothing is also exposed to the risk of contamination.

Model	uvex 3B chem classic	uvex 3B chem classic +	uvex 3B chem light	uvex 4B	uvex 5/6 classic	uvex 5/6 classic light
Art. no.	89880	88498	89843	98375, 89094	98449	17595, 17596
ISO/FDIS 16604/16603 Resistance to contaminated liquids	6 of 6	6 of 6	6 of 6	6 of 6	6 of 6	6 of 6
EN 14126 appendix A Resistance due to mechanical contact with substances containing contaminated liquids	6 of 6	6 of 6	6 of 6	6 of 6	6 of 6	6 of 6
ISO/DIS 22611 Resistance to contaminated liquid aerosols	3 of 3	3 of 3	3 of 3	3 of 3	3 of 3	3 of 3
ISO/DIS 22612 Resistance to contaminated solid particles	3 of 3	3 of 3	3 of 3	3 of 3	3 of 3	3 of 3


Disposable Protection Clothing

Guide to standards and products

Your team's safety at the forefront

From light liquid splashes to hazardous chemicals and radioactive material – employees face various hazards in the course of their daily work. To simplify the process of selecting appropriate protective clothing, the European Union has defined harmonised standards for different protection classes (also called types) for chemical protection – Category III.

The protection type is confirmed by certification and indicates the kind of exposure (dust, liquid, gas) for which the suit is suitable. The attainment of the individual types only ensures the minimum requirements necessary for certification. Please refer to our technical datasheets for details of the individual tests.

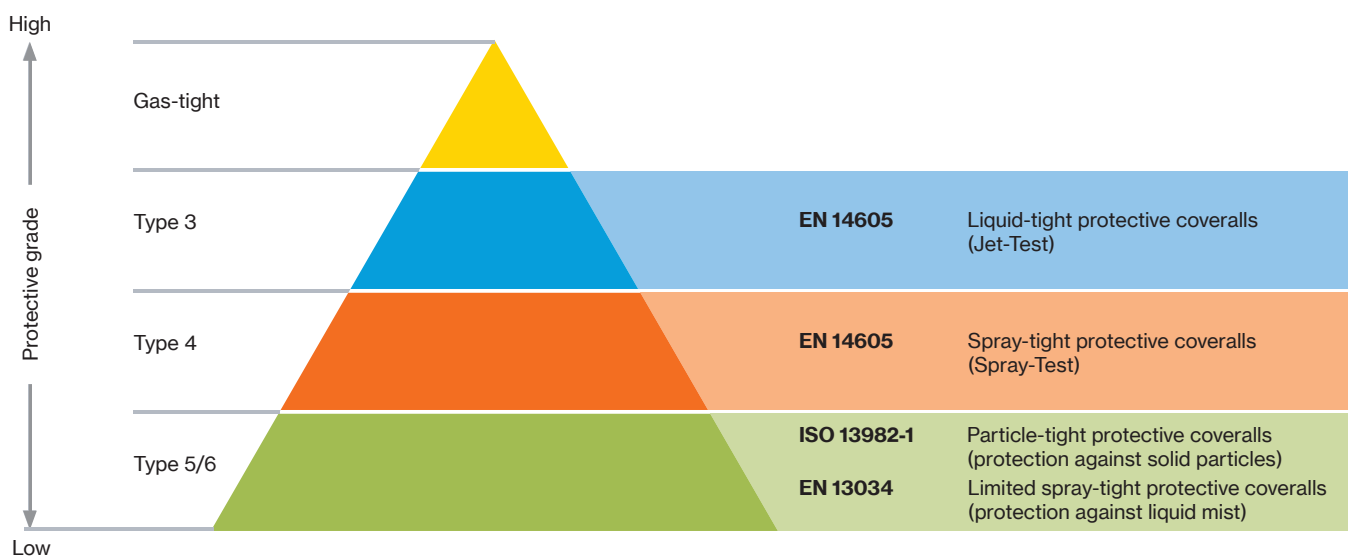

For optimum protection of your team at all times, the breakthrough times of permeation must be taken into account for protective coveralls. uvex protection clothing has undergone the required permeation tests. More information on the breakthrough times of specific chemicals can be found in the technical datasheets or in the uvex Chemical Expert System

Online chemicals database

The uvex Chemical Expert System (CES) offers an extensive chemicals database for choosing the appropriate disposable overall for working with hazardous substances.

As a user, you can create a personal permeation list or receive advice from our specialists. It only takes a few clicks of the mouse to find the right disposable safety clothing and chemical safety gloves to match your specific requirements.



<https://ces.uvex.de>

The breakthrough times were determined under laboratory conditions. As additional stresses, such as higher temperatures and mechanical influences, are often present in practice, this information serves as a guide only. Following contamination, wear or damage, the coverall is to be removed immediately and properly disposed of.

Disposable chemical protection

Guidelines for use

Tips for putting on and taking off



Instructions for putting on and taking off Type 3B (PDF)

Instructions for putting on



Instructions for taking off



Instructions for putting on and taking off Type 5/6 (PDF)

Instructions for putting on/
taking off



Information on selecting the correct size of disposable protective suit

To ensure an optimum fit and keep the wearer safe when working with hazardous substances, uvex disposable overalls are available in multiple sizes. However, the correct size may not be the same as your usual clothing size. Use the following overview to find the correct fit based on your height and chest circumference.

Body height in cm	Chest circumference in cm	Size to choose
164 - 170 cm	84 - 92 cm	S
170 - 176 cm	92 - 100 cm	M
176 - 182 cm	100 - 108 cm	L
182 - 188 cm	108 - 116 cm	XL
188 - 194 cm	116 - 124 cm	XXL
194 - 200 cm	124 - 132 cm	XXXL

Selecting the correct chemical-protection clothing

In order to select appropriate chemical-protection clothing for a specific application within your business, it is important to have your overalls tested by a trained safety officer. Many chemicals can cause serious or irreversible injury to workers if they are not adequately protected. For this reason, chemical-protection clothing should be selected with great care.

Note: It is the responsibility of the user to decide for themselves whether the protective suit used is suitable and correct for the intended use. uvex accepts no responsibility whatsoever in the event of improper use and application of these products.

Proper handling of uvex disposable protective clothing

Storage


uvex disposable protective clothing must be stored in its original packaging, away from light and in a dry environment at temperatures of 15°C–25°C. The products must be disposed of after use in accordance with the respective legal requirements.

Washing disposable protective suits

Disposable overalls are for single use only and must not be washed.

Durability

It is recommended to use the product within five years from the production date, which can be found on the label.

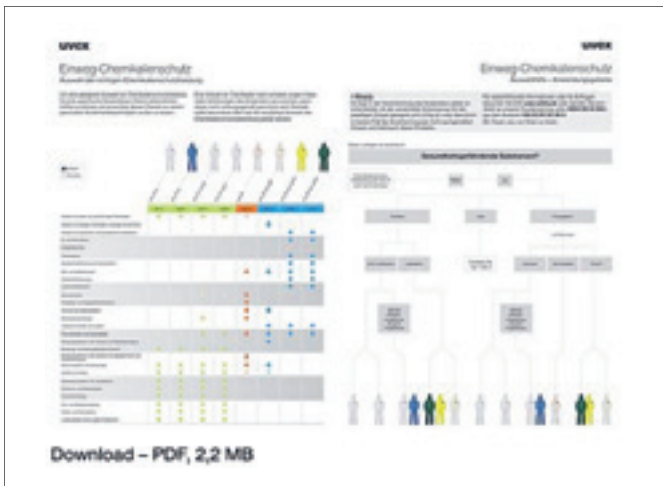

uvex-safety.com/en/overalls