

KIMTECH™

Kimtech™ A5 Sterile Boot Cover with Grasp Ties

Donning Technique

- 1** Select your boot cover size; inspect package for vacuum seal; tear a long notch toward the product to open; flex package to separate boot covers.
- 2** Pull out one boot cover by the cuff.
- 3** Grasp ties with one hand and make sure they don't touch the floor.
- 4** Open up boot cover; point toe toward opening and pull up over calf. Hold ties to prevent touching the floor.
- 5** Hold up foot or rest on bench. Grasp ties at front of boot cover and wrap over shoe/boot; then wrap ties behind ankle and tie in front.
- 6** Place donned boot on the clean side.
- 7** Pull the remaining boot cover out of package by the cuff and repeat steps 3, 4 and 5. Proceed with final gowning steps.



For more information visit us at www.eu.fishersci.com or contact your local Fisher Scientific Representative.

©/™ Trademarks of Kimberly-Clark Worldwide, Inc. ©2020 KCWW Publication code ID 5372.01 EN 06.20

Distributed by Fisher Scientific. Contact us today:

Austria: fishersci.at Belgium: fishersci.be Denmark: fishersci.dk
Germany: fishersci.de Ireland: fishersci.ie Italy: fishersci.it
Finland: fishersci.fi France: fishersci.fr Netherlands: fishersci.nl
Norway: fishersci.no Portugal: fishersci.pt Spain: fishersci.es
Sweden: fishersci.se Switzerland: fishersci.ch UK: fishersci.co.uk

© 2022 Thermo Fisher Scientific Inc. All rights reserved.
Trademarks used are owned as indicated at fishersci.com/trademarks.

 **fisher scientific**
part of Thermo Fisher Scientific