



Kimtech™ A5 Sterile Hood

Donning Technique

- 1** Select your hood; inspect package for vacuum seal; tear along notches toward the product to open.
- 2** Pull out hood and unfold to locate face opening.
- 3** Push one hand through opening while pulling down flaps with other hand.
- 4** Position hood to pull over your head.
- 5** Tilt head up, reach back and pull down flap in back.
- 6** Pull front flap to a comfortable position; if hood has ties, pull back and tie.
- 7** Complete donning with cleanroom mask and goggles.



For more information visit us at www.eu.fishersci.com or contact your local Fisher Scientific Representative.

©/™ Trademarks of Kimberly-Clark Worldwide, Inc. ©2020 KCWW Publication code ID 5372.01 EN 06.20

Distributed by Fisher Scientific. Contact us today:

Austria: fishersci.at Belgium: fishersci.be Denmark: fishersci.dk
Germany: fishersci.de Ireland: fishersci.ie Italy: fishersci.it
Finland: fishersci.fi France: fishersci.fr Netherlands: fishersci.nl
Norway: fishersci.no Portugal: fishersci.pt Spain: fishersci.es
Sweden: fishersci.se Switzerland: fishersci.ch UK: fishersci.co.uk

© 2022 Thermo Fisher Scientific Inc. All rights reserved.
Trademarks used are owned as indicated at fishersci.com/trademarks.